


My Pipette™ Creator App Quick Guide

How to get connected

Thermo Scientific™ E1-ClipTip™ pipettes with software version 1.28 or after


1. Create Thermo Fisher Cloud account

- Go to <http://www.thermofisher.com/cloud> to create your personal Cloud account and sign in.
- Find **My Pipette™ Creator App** from Thermo Fisher Cloud AppConnect homepage and open it

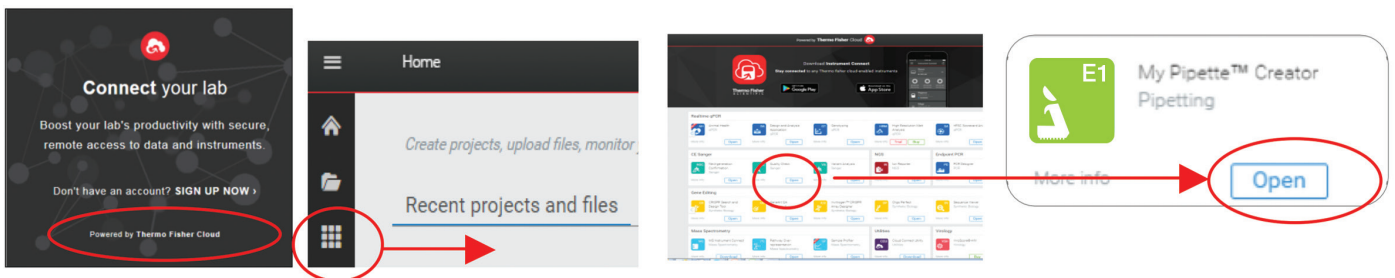

2. Install My Pipette Connect Utility*

- Click 'Add a new E1-ClipTip pipette' function in the My Pipette Creator app.
- Download and install the My Pipette Connect Utility.
- A green icon appears on notification area of PC's taskbar when the utility is installed and running.

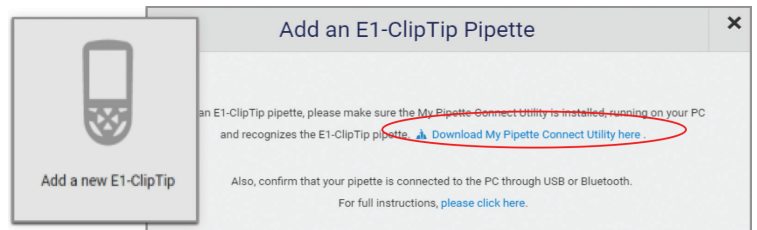


3. Connect E1-ClipTip to PC

- Connect the E1-ClipTip to your PC by:
 - a. Turning on Bluetooth connectivity on your E1-ClipTip pipette** and pairing the pipette to the PC using the Bluetooth Devices function on your PC OR
 - b. Plugging in USB*** cable to the pipette's connector socket.
- A notification pop-up appears on the PC's taskbar and a cloud icon appears on the screen of the E1-ClipTip when the pipette is recognized by the PC.



4. Link your E1-ClipTip to My Pipette Creator app

- Type in the **linking code** provided by the E1-ClipTip pipette (My Pipette >> Thermo Fisher Cloud), and select 'Continue'.
- The pipette will show up under the 'My Pipettes' module on your My Pipette Creator homepage

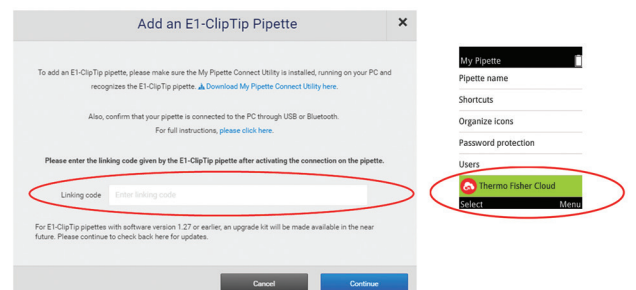

For more information, go to www.thermofisher.com/mypipette

*Installation is required only once on first time usage for specific PC

**only for Bluetooth enabled E1-ClipTip models.

***a USB 2.0 Type-A to USB Micro B plug cable is required

© 2017 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

This product and/or its use in certain methods is/are the subject of patent rights, including US Patent No. 6,778,917, owned by EPPENDORF AG. The product becomes licensed under the aforementioned patents only upon your registration of the product with the Thermo Fisher Cloud.

1517740-01 0217

ThermoFisher
SCIENTIFIC